

AVTECH ELECTROSYSTEMS LTD.

NANOSECOND WAVEFORM ELECTRONICS SINCE 1975

P.O. BOX 265 OGDENSBURG, NY U.S.A. 13669-0265 TEL: 888-670-8729 (USA & Canada) or +1-613-686-6675 (Intl) FAX: 800-561-1970 (USA & Canada) or +1-613-686-6679 (Intl)

BOX 5120, LCD MERIVALE OTTAWA, ONTARIO CANADA K2C 3H4

info@avtechpulse.com - http://www.avtechpulse.com/

INSTRUCTIONS

MODEL AVO-9A-MTB-P1B-T1B-P

200 mA, 200 ps RISE TIME

HIGH PERFORMANCE LASER DIODE DRIVER MODULE

WITH PLUG-IN SOCKET OUTPUT

FOR BUTTERFLY-PACKAGED LASER DIODES

SERIAL	NUMBER:		
		 	$\overline{}$

WARRANTY

Avtech Electrosystems Ltd. warrants products of its manufacture to be free from defects in material and workmanship under conditions of normal use. If, within one year after delivery to the original owner, and after prepaid return by the original owner, this Avtech product is found to be defective, Avtech shall at its option repair or replace said defective item. This warranty does not apply to units which have been dissembled, modified or subjected to conditions exceeding the applicable specifications or ratings. This warranty is the extent of the obligation assumed by Avtech with respect to this product and no other warranty or guarantee is either expressed or implied.

TECHNICAL SUPPORT

Phone: 888-670-8729 (USA & Canada) or +1-613-686-6675 (International) Fax: 800-561-1970 (USA & Canada) or +1-613-686-6679 (International)

E-mail: info@avtechpulse.com World Wide Web: http://www.avtechpulse.com

TABLE OF CONTENTS

WARRANTY	2
TECHNICAL SUPPORT	2
TABLE OF CONTENTS	3
INTRODUCTION	4
SPECIFICATIONS	5
ORIGINAL QUOTATION	6
REGULATORY NOTES	9
FCC PART 18	9
EC DECLARATION OF CONFORMITY	9
DIRECTIVE 2002/95/EC (RoHS)	10
DIRECTIVE 2002/96/EC (WEEE)	10
TERMINAL CONNECTORS	11
SIGNAL CONNECTORS	12
INSTALLING THE DEVICE UNDER TEST	13
NORMAL TEST ARRANGEMENT	15
THERMAL CONTROL (-T1B UNITS)	16
AMPLITUDE CONTROL	16
PROTECTING YOUR INSTRUMENT	17
DO NOT EXCEED 1 MHz	17
PERFORMANCE CHECKSHEET	18

 $Manual\ Reference: \ / fileserver/officefiles/instructword/avo-9/AVO-9A-MTB-P1B-T1B-P, ed 2. odt.$ Last modified November 8, 2010.
Copyright © 2010 Avtech Electrosystems Ltd, All Rights Reserved.

INTRODUCTION

The AVO-9A-MTB-P1B-T1B-P is a high performance DC-powered laser diodr driver module capable of generating up to +200 mA of current into diode loads, at repetition rates up to 1 MHz. The pulse width is adjustable from 0.4 to 4 ns. The rise and fall times are 200 ps or less.

The AVO-9A-MTB-P1B-T1B-P module provides a socket into which the user's laser diode may be inserted. The module internally generates voltage pulses of between 0 and +13V. The module contains the necessary elements to match the laser diode to the pulse generator mainframe. A DC bias current of 0 to +100 mA may be applied to the laser diode by applying the desired DC current to a terminal pin on the module. The module includes an SMA output connector that provides an attenuated coincident replica of the diode current.

The AVO-9A-MTB-P1B-T1B-P must be triggered by an externally-generated TTL-level pulse.

This instrument is intended for use in research, development, test and calibration laboratories by qualified personnel.

SPECIFICATIONS

Model:	AVO-9A-MTB-P1B-T1B-P	
Amplitude, standard:	+20 to +200 mA	
Allowed load voltage range:	0 to 3V	
Pulse width (FWHM):	0.4 - 4 ns	
Duty cycle limits:	N/A	
Minimum PRF:	1 Hz	
Maximum PRF:	1 MHz	
Rise / fall times (20%-80%):	≤ 200 ps	
Related 50Ω series:	AVP-AV-1	
Polarity:	Positive	
Propagation delay:	≤ 150 ns (Ext trig in to pulse out)	
Jitter:	± 35 ps ± 0.015% of sync delay (Ext trig in to pulse out)	
DC offset or bias insertion:	Apply required DC bias current in the range of \pm 100 mA to solder terminal on output module.	
Trigger required: (ext trig mode)	+3 to +5V (TTL), ≥ 50 ns	
Connectors: Out: Other:	-P1B socket, -T1B socket Trig, Monitor: SMA	
Power requirements:	+24V DC and -24V DC	
Dimensions (H×W×D):	43 mm x 76 mm x 200 mm	
Temperature range:	+5°C to +40°C	

ORIGINAL QUOTATION

Date: Tue, 19 Feb 2008 16:28:41 -0500

From: Avtech Sales

To: xxxxx

Subject: Re: Pulse generator

xxxxx,

Following our telephone conversation, I am pleased to quote as follows:

Quote number: 14190

Model number: AVO-9A-MTB-P1B-T1B-P

Description: Laser Diode Driver Module

-MTB special option: The instrument is provided in an DC-powered

externally-triggered module format.

DC power required: +24V and -24V DC

Amplitude: +20 to +200 mA, adjustable

Amplitude control: The amplitude is controlled by a 0 to +10V DC control

voltage applied to one pin of a multi-square-pin header.

Pulse width (FWHM): 0.4 to 4 ns, adjustable.

Pulse width control: The pulse width is controlled by a 0 to +10V DC control voltage applied to one pin of a multi-square-pin header.

Maximum PRF: 1 MHz. The output PRF equals the PRF of the external trigger.

Rise / fall times (20%-80%): < 200 ps

Sync output, Gate input, GPIB/RS-232 ports: none

Monitor output: The standard "MI" output is provided. No "MV" output is provided.

Approximate dimensions (H x W x D): 43 mm x 76 mm x 200 mm

External trigger required: A TTL-level pulse, 50 ns or wider. The input connector is SMA female, and the input impedance is 50 Ohms.

Other: the electrical output is similar to that provided by the standard AVO-9A models described at http://www.avtechpulse.com/laser/avo-9a.

-P1B option: Specifies this common diode configuration (typical of Lumics diodes, and others): The socket will accept pins 8-14 of a standard butterfly package with 0.5 mm wide pins. A positive pulse will be applied to the diode anode (pin 10). Pins 8-9 and 11-13 will be grounded. Four mounting holes on a 8.9 x 26 mm grid will be provided. The diode parasitic resistance (dV/dI at lasing) must be 0 to 2 0hms.

-T1B option: Adds a secondary (low-bandwidth) slide-on socket board for pins 1-7 of a butterfly package. A flexible cable connects the slide-on socket to the output module. A male DB-9 connector is provided on the

output module, which provides access to the thermal control pins of the diode. DB-9 pin 2 connects to diode pin 2 (TH). DB-9 pin 3 connects to diode pin 5 (TH). DB-9 pin 4 connects to diode pin 1 (TEC+). DB-9 pin 5 connects to diode pin 14 (TEC-). The remaining DB-9 pins are unconnected. Pins 3, 4, 6, and 7 of the diode are grounded. Access to the photodiode, if present, is not provided. This option is designed for compatibility with Thorlabs temperature controllers and Lumics laser diodes. It may be suitable for others as well.

Price (Quantity 1-2): \$xxxxx US each, Ex-works Ottawa, Canada.

Price (Quantity 3-4): \$xxxxx US each, Ex-works Ottawa, Canada.

Price (Quantity 5-7): \$xxxxx US each, Ex-works Ottawa, Canada.

Price (Quantity 8-10): \$xxxxx US each, Ex-works Ottawa, Canada.

Quote valid for: 8 weeks

Estimated delivery: 8-12 weeks after receipt of order (excluding export permit* delays).

*Export Permit: These instruments are very high performance pulse generators, which are considered to be "Nuclear-Related Dual-Use Goods" under government regulations. As such, an "End Use Statement" must be completed when ordering. The necessary form is attached (in PDF format). We will use the information in the completed form to apply for an export license from the Canadian government, which will take 1 to 6 weeks to obtain. We cannot ship your order without the license. Please return the completed form to us by fax.

WEEE/ROHS notes: The form available at http://www.avtechpulse.com/green/weee_agreement.pdf must be completed and faxed to Avtech prior to any direct sales by Avtech to European customers.

You will need to confirm that the $\mbox{-P1B}$ and $\mbox{-T1B}$ sockets match your diode pinout.

Please call or email me if I can be of further assistance.

Thank you for your interest in our products!

Regards, Dr. Michael J. Chudobiak Chief Engineer

--- Avtech Electrosystems Ltd. ----- since 1975 ---

PO Box 265 ph: 888-670-8729 or 613-226-5772 Box 5120 Ogdensburg fax: 800-561-1970 or 613-226-2802 LCD Merivale Ottawa, Ontario USA 13669-0265 email: info@avtechpulse.com Canada K2C 3H4 http://www.avtechpulse.com/

Pulse Generators - Laser Diode Drivers - HV Amplifiers Monocycle Generators - Impulse Generators - Pulse Amplifiers Current Pulsers - Function Generators - Frequency Dividers - and more!

REGULATORY NOTES

FCC PART 18

This device complies with part 18 of the FCC rules for non-consumer industrial, scientific and medical (ISM) equipment.

This instrument is enclosed in a rugged metal chassis and uses a filtered power entry module (where applicable). The main output signal is provided on a shielded connector that is intended to be used with shielded coaxial cabling and a shielded load. Under these conditions, the interference potential of this instrument is low.

If interference is observed, check that appropriate well-shielded cabling is used on the output connectors. Contact Avtech (info@avtechpulse.com) for advice if you are unsure of the most appropriate cabling. Also, check that your load is adequately shielded. It may be necessary to enclose the load in a metal enclosure.

If any of the connectors on the instrument are unused, they should be covered with shielded metal "dust caps" to reduce the interference potential.

This instrument does not normally require regular maintenance to minimize interference potential. However, if loose hardware or connectors are noted, they should be tightened. Contact Avtech (info@avtechpulse.com) if you require assistance.

EC DECLARATION OF CONFORMITY



We

Avtech Electrosystems Ltd. P.O. Box 5120, LCD Merivale Ottawa, Ontario Canada K2C 3H4

declare that this pulse generator meets the intent of Directive 89/336/EEC for Electromagnetic Compatibility. Compliance pertains to the following specifications as listed in the official Journal of the European Communities:

EN 50081-1 Emission

EN 50082-1 Immunity

and that this pulse generator meets the intent of the Low Voltage Directive 72/23/EEC as amended by 93/68/EEC. Compliance pertains to the following specifications as listed in the official Journal of the European Communities:

EN 61010-1:2001 Safety requirements for electrical equipment for measurement, control, and laboratory use

DIRECTIVE 2002/95/EC (RoHS)

This instrument is exempt from Directive 2002/95/EC of the European Parliament and of the Council of 27 January 2003 on the Restriction of the use of certain Hazardous Substances (RoHS) in electrical and electronic equipment. Specifically, Avtech instruments are considered "Monitoring and control instruments" (Category 9) as defined in Annex 1A of Directive 2002/96/EC. The Directive 2002/95/EC only applies to Directive 2002/96/EC categories 1-7 and 10, as stated in the "Article 2 - Scope" section of Directive 2002/95/EC.

DIRECTIVE 2002/96/EC (WEEE)


European customers who have purchased this equipment directly from Avtech will have completed a "WEEE Responsibility Agreement" form, accepting responsibility for WEEE compliance (as mandated in Directive 2002/96/EC of the European Union and local laws) on behalf of the customer, as provided for under Article 9 of Directive 2002/96/EC.

Customers who have purchased Avtech equipment through local representatives should consult with the representative to determine who has responsibility for WEEE compliance. Normally, such responsibilities with lie with the representative, unless other arrangements (under Article 9) have been made.

Requirements for WEEE compliance may include registration of products with local governments, reporting of recycling activities to local governments, and financing of recycling activities.



TERMINAL CONNECTORS

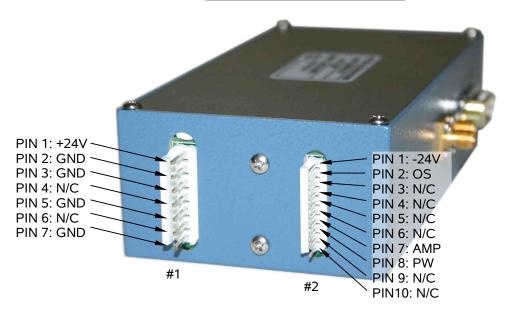

Connectors #1 and #2 are shown above. These supply the DC power and the amplitude and pulse width signals to the module, as well as any required DC offset for the output.

- +24V must be connected to pin 1 of connector #1.
- -24V must be connected to pin 1 of connector #2.

Ground must be connected to pin 2 of connector #1.

The output amplitude is controlled by the voltage (in the range of 0 to +10V) applied to pin 7 of connector #2. The input impedance is several kilohms or higher.


The output pulse width is controlled by the voltage (in the range of 0 to +10V) applied to pin 8 of connector #2. The input impedance is several kilohms or higher.

A forward DC bias may be applied to the laser diode by connecting a DC potential of 0 to +10 Volts to the OS terminal, pin 2 of connector #2. The application of a small forward bias often yields a more ideal diode current waveform (as observed on the MI port).

Connector #1 is an AMP 640445-7 connector. This connector is shipped with a mating AMP 640600-7 insulation-displacement connector installed, for the convenience of the user. The 640600-7 is designed to accept AWG #20 hookup wire.

Connector #2 is an AMP 1-640456-0 connector. This connector is shipped with a mating AMP 4-640621-0 insulation-displacement connector installed, for the convenience of the user. The 4-640621-0 is designed to accept AWG #24 hookup wire.

SIGNAL CONNECTORS

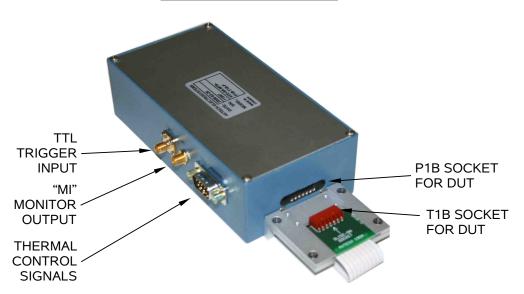

The externally-generated TTL trigger signal must be applied to the SMA connector shown above. The input impedance is 50 Ohms.

An oscilloscope may be used to monitor the MI monitor outputs, as discussed in later sections.


INSTALLING THE DEVICE UNDER TEST

The module will have a "P1B" high-speed socket for pins 8-14 of the diode under test. If the "-T1B" option has been specified, a slide-on socket for pins 1-7 of the diode will also be present.


The "P1B" socket consists of seven high-bandwidth pin sockets. This socket arrangement will accept pins 8-14 of a standard butterfly package with 0.5 mm wide pins. A positive pulse will be applied to the diode anode (pin 10). Pins 8-9 and 11-13 will be grounded.

The optional "T1B" socket consists of a low-bandwidth slide-on socket board for pins 1-7 of a butterfly package. A flexible cable connects the slide-on socket to the output module. A male DB-9 connector is provided on the output module, which provides access to the thermal control pins of the diode. DB-9 pin 2 connects to diode pin 2 (TH). DB-9 pin 3 connects to diode pin 5 (TH). DB-9 pin 4 connects to diode pin 1 (TEC+). DB-9 pin 5 connects to diode pin 14 (TEC-). The remaining DB-9 pins are unconnected. Pins 3, 4, 6, and 7 of the diode are grounded. Access to the photodiode, if present, is not provided. This option is designed for compatibility with Thorlabs temperature controllers and Lumics laser diodes. It may be suitable for others as well.


With no diode installed, the module will look similar to this:

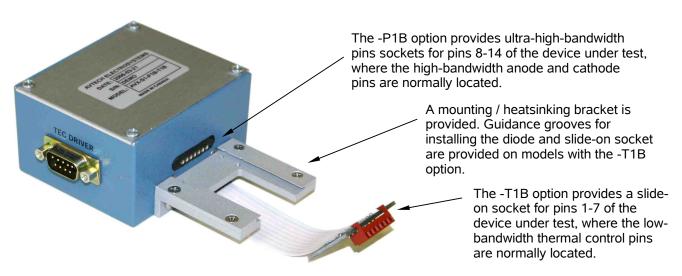

The diode is first installed by sliding pins 8-14 into the "P1B" pin sockets, as shown below:

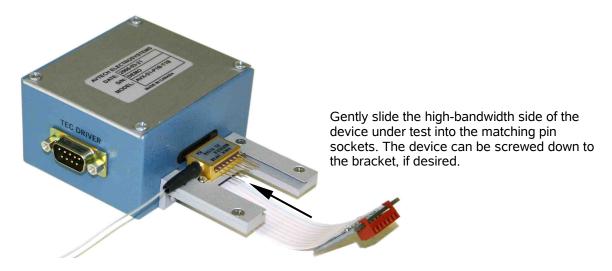


If present, the T1B slide-on socket assembly can then be slid onto pins 1-7, as shown below:

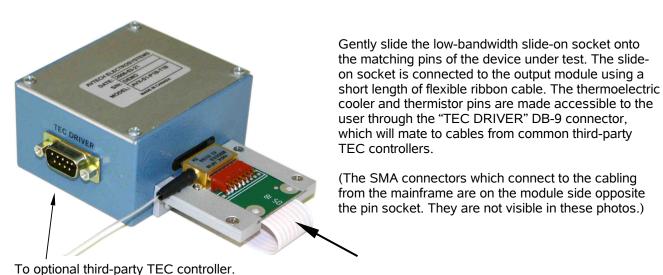

NORMAL TEST ARRANGEMENT

To fully test the instrument, and for normal operation, the module must be connected as shown below. The basic functional equivalent circuit of the output module is shown below:

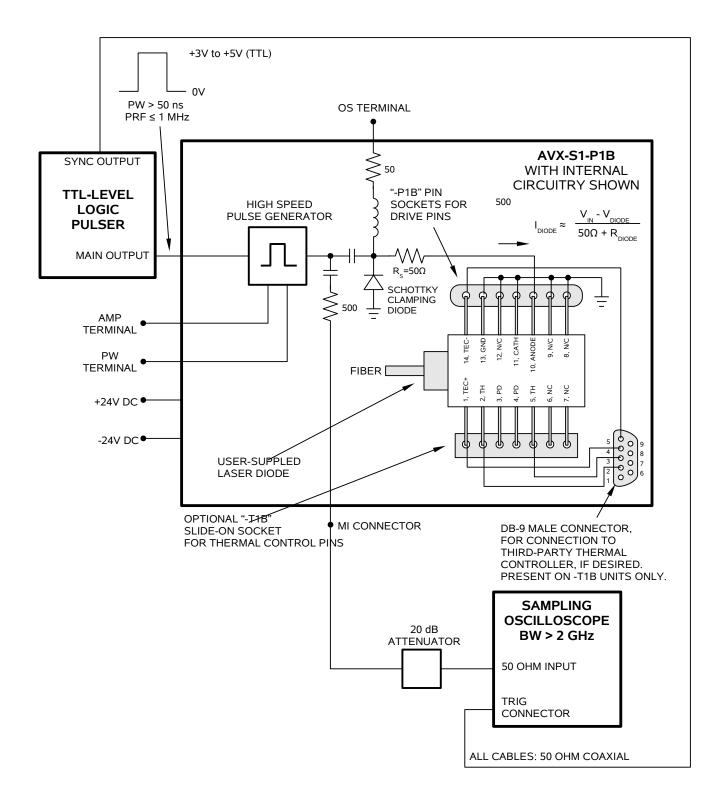

THERMAL CONTROL (-T1B UNITS)

Models with the "-T1B" option include a DB-9 male connector that will mate to third-party temperature controllers. These third-party controller provide a means of controlling the thermoelectric cooler that is typically present in butterfly-packaged laser diodes, if desired. The need for cooling is dependent on the user's application. Cooling is generally recommended by device manufacturers.

AMPLITUDE CONTROL

The pulse current through the diode load installed in the module is given by:

$$I_{DIODE} = (V_{SET} - V_{DIODE}) / (50\Omega + R_{DIODE})$$

where V_{SET} is the amplitude of the internal high-speed voltage pulse generator (between 0 and 13V), V_{DIODE} is the forward voltage drop across the diode (typically 2 or 3V), and R_{DIODE} is typically negligible (compared to the 50 Ohm series resistance).

PROTECTING YOUR INSTRUMENT

DO NOT EXCEED 1 MHz

The output stage may be damaged if triggered by an external signal at a pulse repetition frequency greater than 1 MHz.

PERFORMANCE CHECKSHEET